


Amaxa™ 96-well Shuttle™ Protocol for Mouse ES Cells

Cell Description

This protocol is designed for mouse embryonic stem (ES) cells. Cells derived from mouse blastocysts; round cells, growing in clumps.

Example for Nucleofection™ of Mouse Embryonic Stem Cells

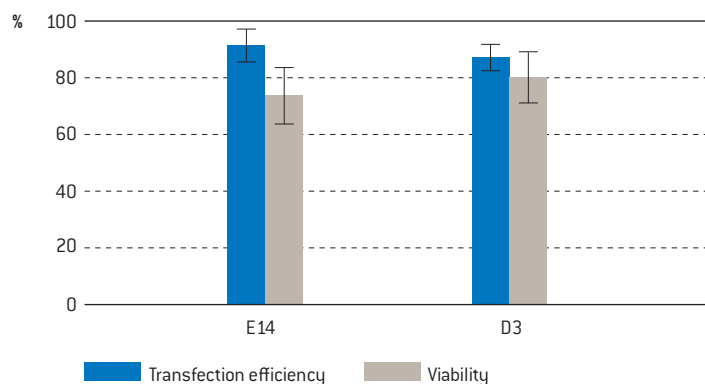

Average transfection efficiency of mouse ES cells 24 hours post Nucleofection™. 5×10^4 mouse ES cells were transfected with program 96-CG-104 and 0.4 μg of pmaxGFP™ Vector. 24 hours post Nucleofection™ cells were analyzed on a FACSCalibur™ with HTS option. Cell viability is given in percent compared to non-transfected control.

Product Description

Recommended Kits

P3 Primary Cell 96-well Nucleofector™ Kits

Cat. No.	V4SP-3096
Size (reactions)	1×96
P3 Primary Cell 96-well Nucleofector™ Solution	2.25 ml
Supplement	0.5 ml
pmaxGFP™ Vector (1 $\mu\text{g}/\mu\text{l}$ in 10 mM Tris pH 8.0)	50 μg
Nucleocuvette™ Plate (s)	1

Cat. No.	V4SP-3960
Size (reactions)	10×96
P3 Primary Cell 96-well Nucleofector™ Solution	22.5 ml
Supplement	5 ml
pmaxGFP™ Vector (1 $\mu\text{g}/\mu\text{l}$ in 10 mM Tris pH 8.0)	50 μg
Nucleocuvette™ Plate (s)	10

Storage and Stability

Store Nucleofector™ Solution, Supplement and pmaxGFP™ Vector at 4°C. For long-term storage, pmaxGFP™ Vector is ideally stored at -20°C. The expiration date is printed on the solution box. Once the Nucleofector™ Supplement is added to the Nucleofector™ Solution, it is stable for three months at 4°C.

Notes

- This protocol is meant to provide an outline for the handling and the Nucleofection™ of mouse ES cells. It has been optimized for mouse ES cell lines ES-E14TG2a (ATCC®-CRL-1821™) and ES-D3 (ATCC®-1934™) but will work with other cells lines as well.
- Experimental results and cell viability may vary depending on the mouse ES cell line used.
- 96-well Nucleofector™ Solutions can only be used with conductive polymer cuvettes, i.e. in the 96-well Shuttle™ Device and in the 4D-Nucleofector™ System. They are not compatible with the Nucleofector™ II/2b Device.

Required Material

Note

Please make sure that the entire supplement is added to the Nucleofector™ Solution.

- Nucleofector™ 96-well Shuttle System (Nucleofector™ Device, version IIS; 96-well Shuttle™ Device; laptop with 96-well Shuttle™ Software)
- Supplemented 96-well Nucleofector™ Solution at room temperature
- Supplied Nucleocuvette™ Plate (s)
- Supplied pmaxGFP™ Vector, stock solution 1 µg/µl

Note

Volume of substrate solution added to each sample should not exceed 10 % of the total reaction volume (2 µl for 20 µl reactions). For positive control using pmaxGFP™ Vector, please dilute the stock solution to reach the appropriate working concentration.

- Substrate of interest, highly purified, preferably by using endotoxin-free kits; A260 : A280 ratio should be at least 1.8
- Nucleocuvette™ compatible tips: epT.I.P.S. (US/CDN: Eppendorf North America, Cat. No. 2491.431, Rest of World: Eppendorf AG, Cat. No. 0030073.266), Matrix TallTips™ (Matrix Technologies Corp., Cat. No. 7281) or LTS Tips (Rainin Instrument, LLC, Cat. No. SR-L10F, SR/SS-L250S, SR/SS-L300S). Before using other types of pipette tips, please ensure they reach the bottom of the Nucleocuvette™ Wells without getting stuck
- 96-well culture plates or culture plates of your choice
- For detaching cells: 0.5 mg/ml Trypsin and 0.2 mg/ml EDTA in PBS and supplemented culture media or PBS/0.5 % BSA
- Mouse ES culture medium: DMEM (ATCC; Cat. No. 30-2002) supplemented with 10 % FCS (ES Cell Qualified) (Invitrogen; Cat. No. 16141-079), 1000 U/ml ESGRO LIF (Chemicon; Cat. No. ESG1107), 0.1 mM 2-Mercaptoethanol and 1 % Pen/Strep (100 µg/ml streptomycin, 100 U/ml penicillin)
- MEF feeder cells: We recommend using irradiated mouse embryonic fibroblasts (MEF) (STO IRR, ATCC® 56-X™) or STO MEF feeder cells (ATCC® CRL-1503™) inactivated by mitomycin-C treatment (see chapter 1)
- For MEF inactivation: Mitomycin-C (Sigma); Ca²⁺- and Mg²⁺-free HBSS
- MEF Culture Medium: DMEM (ATCC 30-2002) + 10 % FCS (ES Cell Qualified) (Invitrogen 16141-079) + 1 % P/S (100 µg/ml streptomycin, 100 U/ml penicillin)
- Prewarm appropriate volume of mouse ES culture medium to 37°C (200 µl per sample)
- Appropriate number of cells (5×10⁴ cells per sample; lower or higher cell numbers may influence transfection results)

1. Pre Nucleofection™

Cell Culture

Note

Mouse ES cells are cultured on irradiated or inactivated mouse embryonic fibroblast (MEF) feeder cells in gelatine-coated plates.

Culture Conditions for MEFs Before Inactivation

- 1.1 Replace medium 2 to 3 times per week
- 1.2 Cells should be passaged 3 times per week with a subcultivation ratio of 1 : 3 to 1 : 10

Inactivation of MEF Feeder Cells

- 1.3 Prepare a mitomycin-C stock solution 1 mg/ml stock, filter sterilized it and store sterile solution at 4°C protected from light
- 1.4 Expand a vial of MEFs as described by ATCC®
- 1.5 Add 1 mg/ml mitomycin-C stock solution to the medium to a final concentration of 10 µg/ml. Return plates to the incubator for 2 to 3 hours
- 1.6 Rinse plates twice with 10 to 15 ml Ca²⁺- and Mg²⁺-free HBSS
- 1.7 Trypsinize as if passaging
- 1.8 Add an equal volume of MEF culture medium without penicillin/streptomycin
- 1.9 Count and freeze cells

Preparation of Gelatin Coated Plates

- 1.10 Add 0.5 grams gelatin to 500 ml endotoxin-free water (0.1 % solution). Gelatin will not be soluble. Do not use glass bottles that have seen detergent. Glass bottles should be cleaned with NaOH when first obtained, and then dedicated to sterile gelatin solution. Do not allow water or gelatin solution to sit unsterilized for any longer than a couple of hours before autoclaving
- 1.11 Autoclave for 30 minutes. Gelatin will solubilize and remain a liquid. Store at room temperature
- 1.12 At least one hour prior to plating of irradiated MEF, coat plate/flask with gelatin solution by placing at least 100 µL per well of gelatin into a 96-well plate (or 3 ml gelatin solution into a 25 cm² flask). Tilt plates/flask in several directions so that liquid covers the entire surface area
- 1.13 Place plates/flasks into a 37°C incubator overnight. Plates/flasks can remain for longer than one day, however they may dry out
- 1.14 Immediately prior to plating of inactivated MEF (inactivated by either radiation or mitomycin-C treatment), aspirate remaining gelatin solution. Either freshly inactivated MEFs or frozen inactivated MEFs can be used
- 1.15 Plate 2×10⁶ inactivated MEF per 25 cm² flask

Cell Culture of Mouse ES Cells

Note

The culture conditions may vary depending on the cell line used. Please refer to more detailed protocols cited under Additional Information before starting experiments.

- 1.16 The medium should be replaced every day!
- 1.17 Mouse ES cells should be passaged every second day and seeded on on freshly plated feeder cells. We recommend a subcultivation ratio of 1 : 4 to 1 : 10 (1×10^6 mouse ES cells per 25 cm^2 flask)

Trypsinization

- 1.18 24 hours before passaging the mouse ES cells plate inactivated MEF feeder cells on a gelatine-coated flask (2×10^6 inactivated MEF per 25 cm^2 flask)
- 1.19 Detach the mouse ES cells from the plate by trypsinization
- 1.20 Stop trypsinization by adding 4-fold media
- 1.21 As MEF feeder cells are also detached by this treatment, purge the cell suspension into an untreated flask (cell-culture treated) for 1 hour. MEF feeder cells accompanying the mouse ES cells will attach to the surface of the flask during this time
- 1.22 Collect the supernatant containing the mouse ES cells and plate it into a gelatine coated flask with inactivated feeder cells

2. Nucleofection™

Note

Prepare a 96-well culture plate with feeder cells for post Nucleofection™ culturing of the transfected cells 24 hours before Nucleofection™. Therefore plate 6.3×10^4 inactivated feeder cells into each well of a gelatin coated 96-well plate (6×10^6 cells for a complete 96-well plate).

One Nucleofection™ Sample Contains

- 5×10^4 cells
 - 0.4–0.8 µg plasmid DNA (in 1–2 µl H₂O or TE) or 0.4 µg pmaxGFP™ Vector or 30–300 nM siRNA (0.6–6 pmol/sample)
 - 20 µl P3 Primary Cell 96-well Nucleofector™ Solution
- 2.1 Please make sure that the entire supplement is added to the Nucleofector™ Solution
 - 2.2 Start Nucleofector™ 96-well Shuttle™ Software, verify device connection and upload experimental parameter file (for details see device and software manuals)
 - 2.3 Select the appropriate 96-well Nucleofector™ Program **96-CG-104**
 - 2.4 Prepare cell gelatin coated culture plates by filling appropriate number of wells with desired volume of recommended culture media, e.g. 120 µl* (see note at the end of this chapter) for one well of a 96-well plate precoated with gelatin and inactivated feeder and pre-incubate/equilibrate plates in a humidified 37°C/5% CO₂ incubator

- 2.5 Pre-warm an aliquot of culture medium to 37°C (80 µl per sample*)
- 2.6 Prepare 0.4–0.8 µg plasmid DNA or 0.4 µg pmaxGFP™ Vector or 30–300 nM siRNA (0.6–6 pmol/sample)
- 2.7 Prepare the required amount of mouse ES cells by trypsinization as described in 1.18–1.22
- 2.8 After incubating the trypsinized cell suspension for 1 hour on uncoated cell culture flasks, collect the mouse ES cells from by centrifugation (125×g, 10 minutes, RT). Remove supernatant completely
- 2.9 Resuspend the cell pellet carefully in 20 µl room temperature 96-well Nucleofector™ Solution per sample

A: One or several substrates (DNAs or RNAs) in multiples

- Prepare mastermixes by dividing cell suspension according to number of substrates
- Add required amount of substrates to each aliquot (max. 2 µl per sample)
- Transfer 20 µl of mastermixes into the wells of the 96-well Nucleocuvette™ Modules

B: Multiple substrates (e.g. Library Transfection)

- Pipette 20 µl of cell suspension into each well of a sterile U- or V-bottom 96-well microtiter plate
- Add 2 µl substrates (maximum) to each well
- Transfer 20 µl of cells with substrates into the wells of the 96-well Nucleocuvette™ Modules

Note

It is advisable to pre-dispense each cell suspension into a sterile round-bottom 96-well plate or to pipet from a pipetting reservoir for multi-channel pipettes. Use a multi-channel or single-channel pipette with suitable pipette tips. As leaving cells in 96-well Nucleofector™ Solution for extended periods of time may lead to reduced transfection efficiency and viability it is important to work as quickly as possible. Avoid air bubbles while pipetting.

- 2.10 Gently tap the Nucleocuvette™ Plate to make sure the sample covers the bottom of the well
- 2.11 Place 96-well Nucleocuvette™ Plate with closed lid into the retainer of the 96-well Shuttle. Well "A1" must be in upper left position
- 2.12 Start 96-well Nucleofection™ Process by either pressing "Upload and start" in the 96-well Shuttle™ Software or pressing "Upload" in the 96-well Shuttle™ Software and then the "Start" button at the 96-well Shuttle™ (for both options please refer to the respective Manual)
- 2.13 After run completion, open retainer and carefully remove the 96-well Nucleocuvette™ Plate from the retainer
- 2.14 Resuspend cells with 80 µl* (recommendation for 96-well plates) or desired volume of pre-warmed medium (maximum cuvette volume 200 µl). Mix cells by gently pipetting up and down two to three times

- 2.15 Plate desired amount of cells in culture system of your choice. Recommendation for 96-well plates: Transfer 80 µl of resuspended cells to 120 µl pre-warmed medium prepared in 96-well culture plates*

*** Note**

The indicated cell numbers and volumes have been found to produce optimal 96-well Nucleofection™ Results in most cases. However, depending on your specific needs you may wish to test an extended range of cell numbers. Cell numbers and volumes can be adapted such that fewer cells are transferred or duplicate plates can be seeded.

3. Post Nucleofection™

- 3.1 Incubate the cells in humidified 37°C/5% CO₂ incubator until analysis. Gene expression or down regulation, respectively, is often detectable after only 8–16 hours

Additional Information

Up-To-Date List of all Nucleofector™ References

www.lonza.com/nucleofection-citations

Technical Assistance and Scientific Support

USA/Canada

Tel 800 521 0390 (toll-free)

Fax 301 845 8338

scientific.support@lonza.com

Europe and Rest of World

Tel +49 221 99199 400

Fax +49 221 99199 499

scientific.support.eu@lonza.com

References

1. Hogan, Constantini and Lacy, Manipulation of the mouse embryo, Cold Spring Harbor laboratory Press
2. David A. Conner Harvard Medical School, Boston, Massachusetts
3. Current Protocols in Molecular Biology Copyright © 2003 John Wiley & Sons, Inc.
4. Frederick M. Ausubel, Roger Brent, Robert E. Kingston, David D. Moore, J.G. Seidman, John A. Smith, and Kevin Struhl (eds.)
5. DOI: 10.1002/0471142727.mb2302s51 Online Posting Date: May, 2001 Print Publication Date: July, 2000

www.lonza.com

Lonza Cologne GmbH – 50829 Cologne, Germany

Please note that the Amaxa™ Nucleofector™ Technology is not intended to be used for diagnostic purposes or for testing or treatment in humans. The Nucleofector™ Technology, comprising Nucleofection™ Process, Nucleofector™ Device, Nucleofector™ Solutions, Nucleofector™ 96-well Shuttle™ System and 96-well Nucleocuvette™ Plates and Modules is covered by patent and/or patent-pending rights owned by Lonza Cologne GmbH. Amaxa, Nucleofector, Nucleofection, 96-well Shuttle, Nucleocuvette and maxGFP are either registered trademarks or trademarks of the Lonza Cologne GmbH in Germany and/or U.S. and/or other countries. FACSCalibur is a trademark of Becton Dickinson. TallTips are a registered trademark of Matrix Technologies Corporation. Other product and company names mentioned herein are the trademarks of their respective owners.

This kit contains a proprietary nucleic acid coding for a proprietary copepod fluorescent protein intended to be used as a positive control with this Lonza product only. Any use of the proprietary nucleic acid or protein other than as a positive control with this Lonza product is strictly prohibited. USE IN ANY OTHER APPLICATION REQUIRES A LICENSE FROM EVROGEN. To obtain such a license, please contact Evrogen at license@evrogen.com. The CMV promoter is covered under U.S. Patents 5,168,062 and 5,385,839 and its use is permitted for research purposes only. Any other use of the CMV promoter requires a license from the University of Iowa Research Foundation, 214 Technology Innovation Center, Iowa City, IA 52242. The use of this product in conjunction with materials or methods of third parties may require a license by a third party. User shall be fully responsible for determining whether and from which third party it requires such license and for the obtainment of such license. No statement is intended or should be construed as a recommendation to infringe any existing patent.

© Copyright 2009, Lonza Cologne GmbH. All rights reserved – D4SP-3011 2011-01